

	SOFTWARE RELEASE NOTE	1(4)
Document:	Software Release Note TPLS 3.3.0	Document: TPLS-3_3_0
Project:	TPLS Software Suite	Revision: 1

1 Revision History

Revision	Date	Implemented by	Details
1	January 2024	Doug Lexvold	Initial Release

2 Review & Approval:

Signature: Mike Dingle, Senior Product Specialist 	Date Jan 10 2024
Signature 	Date Jan 8 2024


	SOFTWARE RELEASE NOTE	2(4)
Document:	Software Release Note TPLS 3.3.0	Document: TPLS-3_3_0
Project:	TPLS Software Suite	Revision: 1

Table of Contents

1 Revision History 1

2 Review & Approval: 1

3 Introduction..... 3

4 Change Detail..... 3

 4.1 New Features..... 3


 4.2 Issues Resolved 3

5 Testing..... 3

6 Compatibility Issues..... 4

7 Installation..... 4

8 Usage Notes 4

	SOFTWARE RELEASE NOTE	3(4)
Document:	Software Release Note TPLS 3.3.0	Document: TPLS-3_3_0
Project:	TPLS Software Suite	Revision: 1

3 Introduction

The purpose of this document is to document the release of software for production of TrakPro Lite Secure (TPLS) v3.3.0. It describes the changes to the software beginning with release 3.2.0.

This release of TPLS is issued to be installed subsequent to the installation of previous versions of TPLS 3.x and as a new version to be installed on systems that have not previously had TPLS installed.

TPLS 3.3.0 has been tested on Windows 10, Windows 11 and Windows Server 2019. Supported operating systems therefore include, Windows 10 (64 bit), Windows 11 (64 bit) and Windows Server 2019.

4 Change Detail

TrakPro Lite Secure 3.3.0 incorporates additional features not available in previous versions of the software.

4.1 New Features

Feature Name	Description
N/A	N/A


4.2 Issues Resolved

TrakPro Lite Secure v3.3.0 is an update to release 3.2.0 and resolves the following issues:

Issue Name	Description
Zone deletion not logged in the audit log.	If a zone was deleted, this action was not being included in the audit log. This action is now logged in the audit log.
Audit Log End Date adds a day	When printing the Audit Log a day was added to the End Date. A day is no longer added to the End Date when printing the Audit Log.
Deleted recipes not included in audit log entries.	The name of a deleted recipe would not get included in the Audit log entry. Now, when a recipe has been deleted, its name will appear in audit log entries where appropriate.
Inconsistent date and time stamps	The date and time formats used in the headers and footers of reports were inconsistent. All date and time stamps are now consistently formatted as dd/mm/yyyy and 24-hour clock.
Prerequisite software is no longer supported.	Some software that was required by the installer is no longer supported. End of Life prerequisites have been removed from the installer.

5 Testing

The changes detailed above have been tested on Windows 10, Window 11 and Windows Server 2019.

	SOFTWARE RELEASE NOTE	4(4)
Document:	Software Release Note TPLS 3.3.0	Document: TPLS-3_3_0
Project:	TPLS Software Suite	Revision: 1

6 Compatibility Issues

This update is compatible with all versions of TPLS currently released and CCOPC instruments running ModbusMap2 version 2.00 (no zone support) and 2.01 (zone support). It also supports the AeroTrak 9303.

7 Installation

Run TrakProLiteSecureSetup-3.3.exe. Full instructions will be given during the installation.

8 Usage Notes

None.